

Physical Specifications

Table 3 Physical Specifications

Type	Specification	Comments
Weight	11 kg (25 lbs)	
Dimensions (width × depth × height)	140 x 345 x 435 mm (5.5 x 13.5 x 17 inches)	
Line voltage	100 – 240 VAC, ± 10%	Wide-ranging capability
Line frequency	50 or 60 Hz, ± 5%	
Power consumption	180 VA, 55 W / 188 BTU	Maximum
Ambient operating temperature	4–55 °C (41–131 °F)	
Ambient non-operating temperature	40–70 °C (-4–158 °F)	
Humidity	< 95%, at 25–40 °C (77–104 °F)	Non-condensing
Operating Altitude	Up to 2000 m (6500 ft)	
Non-operating altitude	Up to 4600 m (14950 ft)	For storing the module
Safety standards: IEC, CSA, UL	Installation Category II, Pollution Degree 2	For indoor use only. Research Use Only. Not for use in Diagnostic Procedures.

Performance Specifications

Table 4 Performance Specification Agilent 1200 Series Quaternary Pump

Type	Specification
Hydraulic system	Dual plunger in series pump with proprietary servo-controlled variable stroke drive, floating plungers and active inlet valve
Setable flow range	0.001 – 10 ml/min, in 0.001 ml/min increments
Flow range	0.2 – 10.0 ml/min
Flow precision	£ 0.07% RSD, or £ 0.02 min SD whatever is greater, based on retention time at constant room temperature
Flow accuracy	± 1% or 10 µl/min whatever is greater
Pressure	Operating range 0 – 40 MPa (0 – 400 bar, 0 – 5880 psi) up to 5 ml/min Operating range 0 – 20 MPa (0 – 200 bar, 0 – 2950 psi) up to 10 ml/min
Pressure pulsation	< 2 % amplitude (typically < 1 %), at 1 ml/min isopropanol, at all pressures > 1 MPa (10 bar)
Compressibility compensation	User-selectable, based on mobile phase compressibility
Recommended pH range	1.0 – 12.5, solvents with pH < 2.3 should not contain acids which attack stainless steel
Gradient formation	Low pressure quaternary mixing/gradient capability using proprietary high-speed proportioning valve Delay volume 800 – 1100 µl, dependent on back pressure
Composition Range	0 – 95 % or 5 – 100 %, user selectable
Composition Precision	< 0.2 % RSD, at 0.2 and 1 ml/min
Control and data evaluation	Agilent control software (e.g. ChemStation, EZChrom, OL, etc.)
Communications	Controller-area network (CAN), GPIB, RS-232C, APG Remote: ready, start, stop and shut-down signals, LAN optional

Table 4 Performance Specification Agilent 1200 Series Quaternary Pump

Safety and maintenance	Extensive diagnostics, error detection and display (through control module and Agilent Lab Monitor & Diagnostic Software), leak detection, safe leak handling, leak output signal for shutdown of pumping system. Low voltages in major maintenance areas.
GLP features	Early maintenance feedback (EMF) for continuous tracking of instrument usage in terms of seal wear and volume of pumped mobile phase with user-settable limits and feedback messages. Electronic records of maintenance and errors.
Housing	All materials recyclable.